



Quality. Cooled. Cameras.



ProLine LDR Radiography Imaging Systems

## **Welcome**

Thank you for purchasing an FLI camera. We know that your new camera will bring you years of enjoyment and excellent imaging results.

This User Guide is intended as a reference tool for you to use with the ProLine LDR Imaging System. Please read it and follow the procedures to ensure trouble-free installation of your hardware and software.

If you have any questions about your purchase, please contact us.

## **Contact Information**

**Finger Lakes Instrumentation, LLC**  
**7287 West Main Street**  
**Lima, New York 14485**

**Web: [www.flicamera.com](http://www.flicamera.com)**  
**Email: [support@flicamera.com](mailto:support@flicamera.com)**  
**Fax: 585-624-9879**  
**Phone: 585-624-3760**

# Contents

ProLine LDR Radiography Imaging Systems .....	1
Welcome.....	2
Contact Information .....	2
Check your Shipment .....	4
Product Safety.....	4
ProLine LDR Imaging System Connections .....	5
Setting-up and Powering Your System .....	6
Appendix A - External Triggering .....	9
Exposure Active Indicator .....	9
Shutter Open Indicator.....	10
Appendix B – Troubleshooting.....	11
Appendix C - Camera Specifications.....	12
ProLine PL09000LDR LINOS System .....	12
ProLine PL16803LDR-LINOS System .....	13
Appendix D - Quantum Efficiency Curves .....	14
Kodak KAF-09000.....	14
Kodak KAF-16803.....	14
Appendix E – Warranty for ProLine Imaging Systems .....	15
Appendix F – Camera and Supplies Return Procedure .....	16
Return Addresses by Carrier .....	17
UPS and Fed Ex Returns.....	17
USPS Returns.....	17

## Check your Shipment

Please ensure that all the components have arrived safely and verify that all the items you ordered were properly delivered to you. In the unlikely event of a missing or damaged component, immediately notify your FLI dealer or FLI.

A camera order should include the following items:

- Camera with shutter cover installed (save this cover)
- Camera Test Report
- USB cable
- 12 volt power supply
- Power supply line cord
- Packing list
- User Guide


If you ordered accessories or options, these items should be included. Please check that your order is complete.

## Product Safety

The ProLine LDR Imaging System is shipped with a 12-volt power supply. Do not use any other power supply with this System or use the power supply in a way other than described in this Guide as it may cause damage to the Camera that will not be covered under the warranty.

If you are concerned about lightning strikes in the area in which you use your Camera, you may want to take safety precautions as electrical surges can damage electrical equipment. We recommend that when your Camera is not in use that you unplug the Camera from power and unplug the USB cable from the Camera.

# ProLine LDR Imaging System Connections

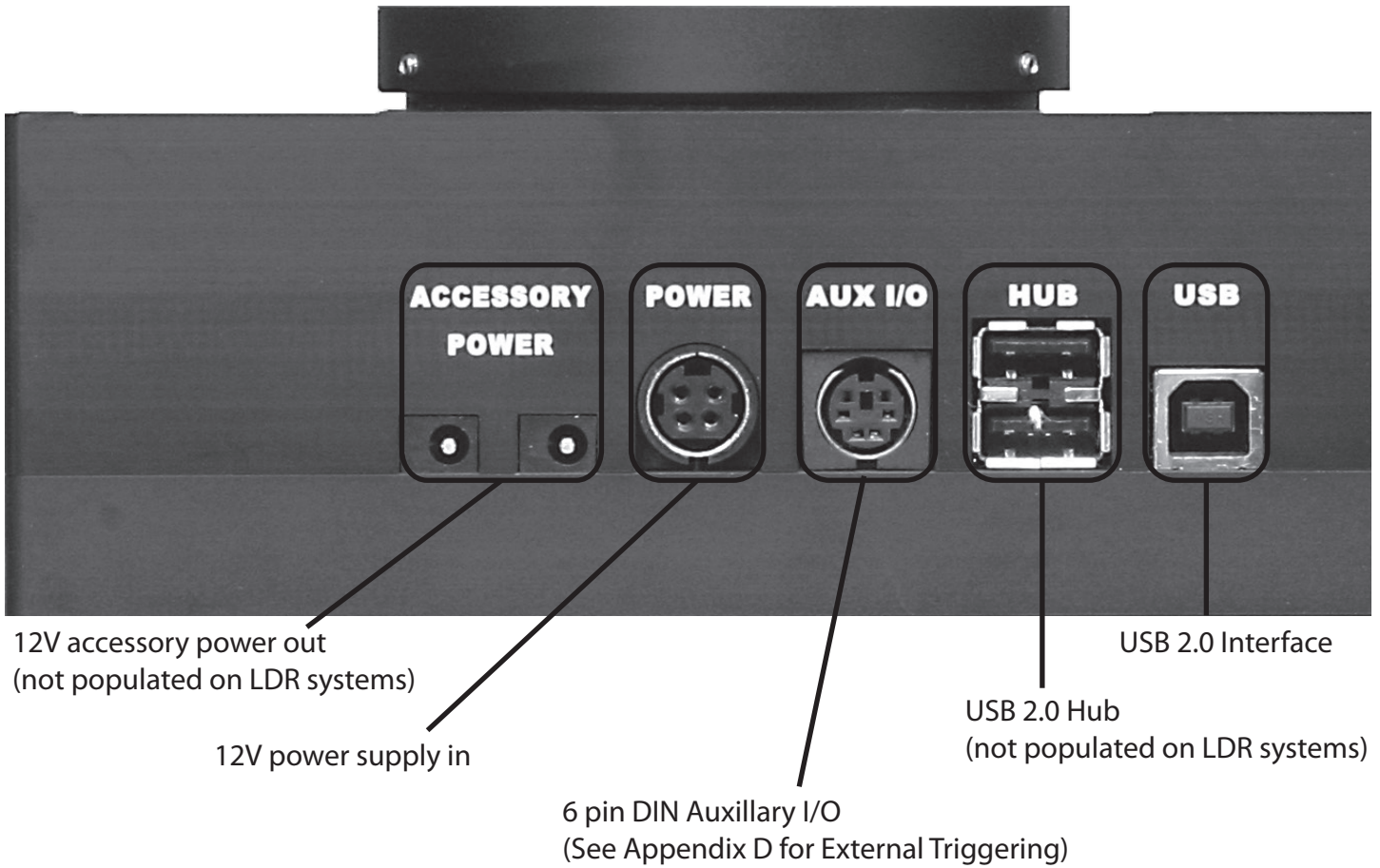

# Setting-up and Powering Your System

Assemble all System components in one area before beginning this set up procedure. The image below depicts an assembled System.


Note: Run the FLI Software Installation Kit before you set up and power your system to ensure the system recognizes and correctly initialized the camera.



Setup the components as described in the procedure below.

1. Attach the USB cable between your ProLine camera and your PC.
2. Connect the 12-volt power supply between the camera's 4-pin "DC Power Jack" connector and a wall socket or switched AC power strip. The camera's fans will start running. Your computer should display a screen indicating that the system has found new hardware.
3. To connect a focuser or a color filter wheel to the camera you will need a short cable. A short cable kit is available from FLI. Install the power cable from the camera to the focuser or from the camera to the color filter wheel. If both are accessories were ordered, two sets of cables are provided.

Note: Refer to Appendix A for information about external triggering using the 6 pin DIN Auxiliary I/O connector.

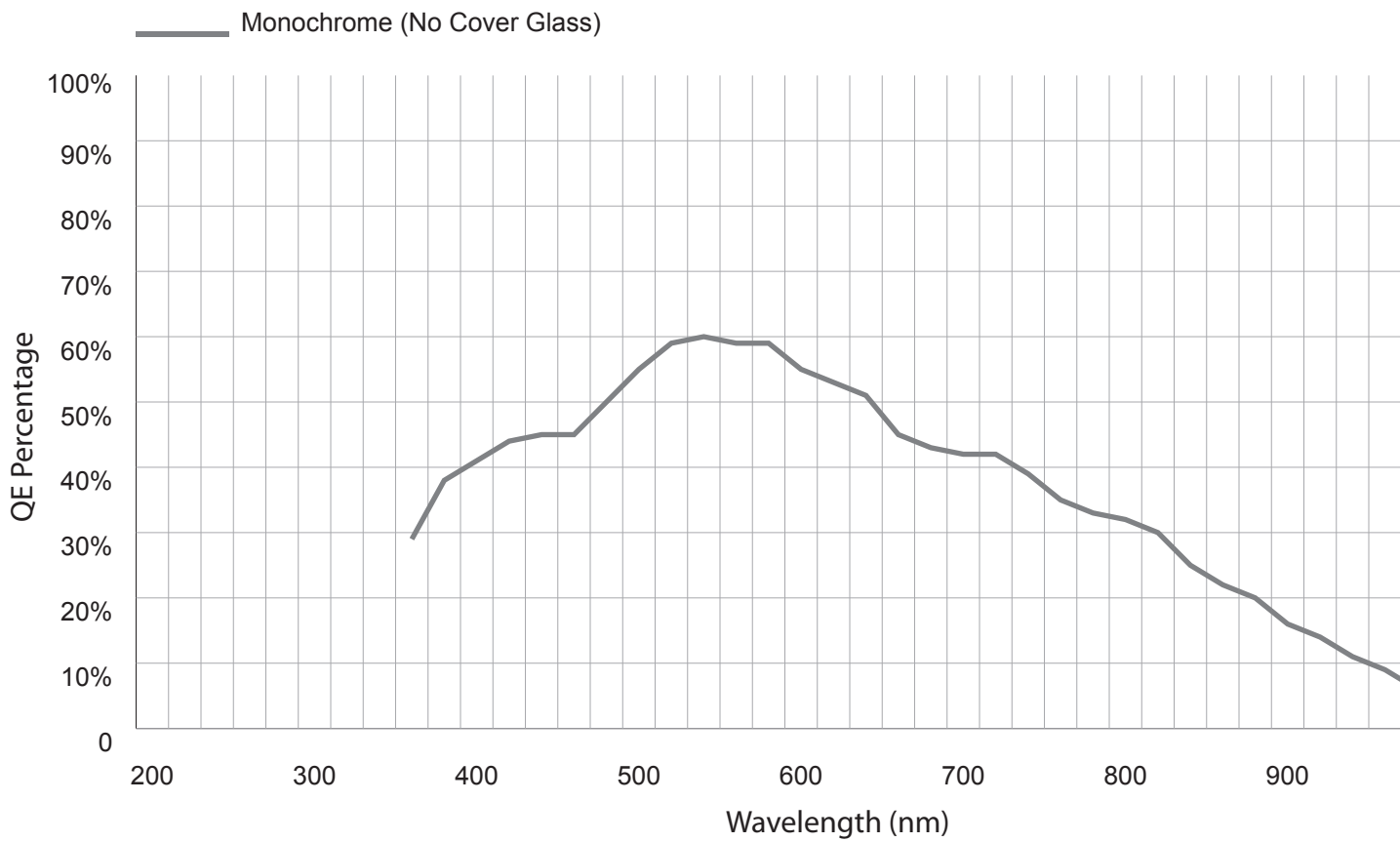


## Camera, Spacer, Lens Configuration

Several cameras, CCD sensors and lenses are available to fit your custom application. A camera unit consists of a camera body with a CCD sensor, a Linos lens and, if necessary, a black aluminum spacer(s). Figure 3 shows the relationship of these components and Table 1 provides information on the specific configurations. In summary:

- The 3801 series camera body can contain one of two different CCD sensors and can accommodate two different types of lenses. Spacers are not required with any 3801 camera body-lens configuration.
- The two PL series camera bodies contain different CCD sensors, yet each can accommodate the three available lenses. Depending on the lens, you will either need no spacer, or you will need one or three spacers stacked on top of each other to ensure the proper back focus distance. If you order camera body-lens unit that requires a spacer(s), it (they) will be installed in the camera body prior to shipping. If you subsequently order a 3801-382 lens, you will be shipped two additional spacers that are required.


### Camera/Lens Spacers

Camera	CCD	# of Spacers	Lens
3801-453-20	9000	0	3801-437-000-20-0102c
3801-453-21	16803	0	3801-437-000-20-0102c
3801-453-22	9000	0	3801-438-000-20-0001a
3801-453-23	16803	0	3801-438-000-20-0001a
PL09000	9000	3	3801-382-000-21_az_d
PL09000	9000	0	3801-437-000-20-0102c
PL09000	9000	1	3801-438-000-20-0001a
PL16803	16803	3	3801-382-000-21_az_d
PL16803	16803	0	3801-437-000-20-0102c
PL16803	16803	1	3801-438-000-20-0001a

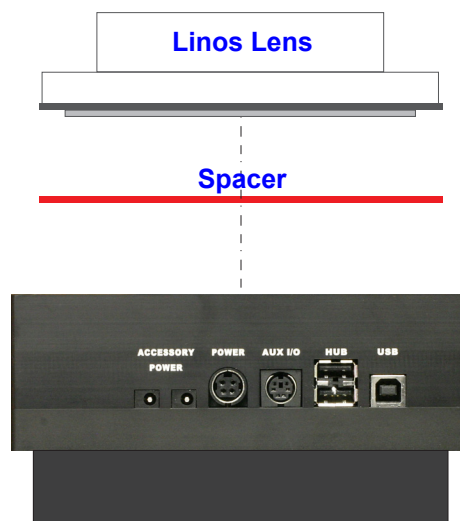

## Appendix A - External Triggering

The ProLine Imaging Systems has a six pin connector shown in the image below.

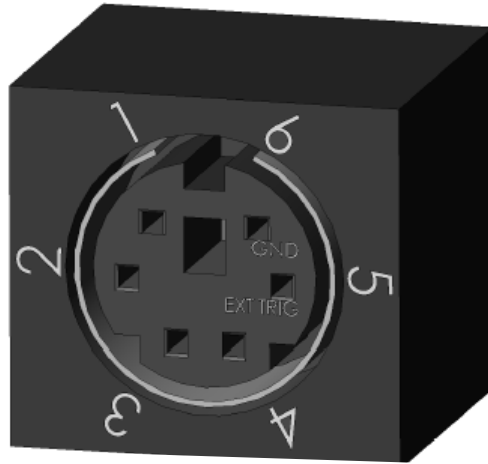

Pin 6 is the electrical ground. Pin 5 is the logic input for the external trigger. (This input has a 22.1K pullup resistor to +3.3V). The external trigger can be setup for either rising edge or falling edge trigger.

To externally trigger the camera using FLIGrab

1. Select the Shutter Control.
2. Select the desired triggering on the pull down menu:
  - ∞ Select “External trigger on LOW” for low edge triggering.
  - ∞ Select “External trigger on HIGH” for high edge triggering.
3. Select “Done”.
4. Open the Grab Menu.
5. Select desired frame and exposure settings.
6. Click the Grab button.
7. The camera will now wait for an external trigger.

A camera waiting for an external trigger “HIGH” will initiate a frame grab when a 0V transition to +3.3V signal or equivalent LVTTTL signal is applied to pin 5.

A camera waiting for an external trigger “LOW” will initiate a frame grab when a +3.3V transition to 0V signal or equivalent LVTTTL signal is applied to pin 5.

### Exposure Active Indicator

---

Pin 1 is an exposure active indicator. It becomes an LVTTTL high with respect to the ground (pin6) when the camera exposure starts and it becomes an LVTTTL low when the camera exposure is finished. It will remain low until the next exposure. It has a 332 Ohm source impedance.

## Shutter Open Indicator

---

Pin 2 on the AUX I/O connector provides an LVTTTL compatible signal for shutter open indication. When the shutter is opened, pin 2 is pulled to +3.3V through an internal 22.1Kohm resistor. When the shutter is closed, pin 2 is pulled down to ground through a 332-ohm resistor within the camera. Pin 6 is the ground reference for this signal.

Pins 3 and 4 are uncommitted camera I/O signals.

### Notes:

The mating part for the external trigger connection is a 6-pin male (CUI MD-60).

The Digi-Key part number for this part is CP-2060-ND.

## Appendix B – Troubleshooting

Issue	Remedy
The fan on the camera does not turn.	Check that the camera is plugged into a suitable power supply and that power is available.
The fan turns but when I use FLIGrab, but the camera does not work or works erratically.	Make certain that all cables are well seated and are not stretched.
FLIGrab runs, the camera fan turns but FLIGrab cannot locate the camera.	Please turn off your computer and camera and leave them switched off for two to three minutes, this will flash the USB chips. Try again.
I get a Windows "Communication" error.	Please turn off your computer and camera and leave them switched off for two to three minutes, this will flash the USB chips. Try again.
Everything is connected and FLIGrab has been installed but the software cannot locate the camera.	<p>Please make certain that you have installed the software and drivers as per the instructions in this manual. Letting Windows locate the camera during installation is likely to create problems.</p> <p>Uninstall FLIGrab and install it again following the installation process.</p>
FLIGrab runs and has identified my camera. I have Grabbed an image but the image appears as a flat black frame.	You are over exposing or under exposing. Check the settings in the software.

# Appendix C - Camera Specifications

## ProLine PL09000LDR LINOS System


<b>CCD Specifications</b>	
Sensor	Kodak KAF-09000
Grade	Standard
CCD Format	3056 (H) x 3056 (V) = 9.3 Mp
Chip Size	38.6 mm (H) x 37.76 mm (V)
Pixel Size	12 $\mu$ m (H) x 12 $\mu$ m (V)
Saturation Signal	120 K e-
Read Noise	7 electrons @ 4 MHz
Dark Current Doubling Temp	7 deg C
<b>Camera Electrical Performance</b>	
Fast Readout Speed	8 MHz
Slow Readout Speed	1 MHz
Typical Readout Noise @ 8 MHz	12-14 electrons RMS
Typical Readout Noise @ 1 MHz	8-10 electrons RMS
Typical Cooling Performance	50 deg C below ambient temperature
I/O Control	External Trigger Control is LVTTTL
Non-Linearity	<1% of signal
Anti-dew	Standard
<b>Camera Physical Characteristics</b>	
Weight	5.5 lbs
HWD	4" x 6.2" x 6.2"
Sealed Chamber	CCD
Power Requirements	12V
Window	1 mm quartz
Back Focus Distance (CCD to Lens)	6.59mm
Mounting Method	4-40 UNC
<b>Lens: Ultimate Imaging Quality under Extreme Operating Conditions</b>	
Aperture	1:1.0 in (Lowest Possible Exposure Times)
Objective Field	43cm x 43cm

## ProLine PL16803LDR-LINOS System


<b>CCD Specifications</b>	
Sensor	Kodak KAF-16803
Grade	Standard
CCD Format	4096 (H) x 4096 (V) = 16.8 Mp
Chip Size	38.6 mm (H) x 37.76 mm (V)
Pixel Size	9 $\mu$ m (H) x 9 $\mu$ m (V)
Saturation Signal	85 K e-
Read Noise	9 electrons @ 4 MHz
Dark Current Doubling Temp	6.6 deg C
<b>Camera Electrical Performance</b>	
Fast Readout Speed	8 MHz
Slow Readout Speed	1 MHz
Typical Readout Noise @ 8 MHz	12-14 electrons RMS
Typical Readout Noise @ 1 MHz	8-10 electrons RMS
Typical Cooling Performance	Up to 45-50 deg C below ambient temperature
I/O Control	External Trigger Control is LVTTTL
Non-Linearity	<1% of signal
Anti-dew	Standard
<b>Camera Physical Characteristics</b>	
Weight	5.5 lbs
HWD	4" x 6.2" x 6.2"
Sealed Chamber	CCD
Power Requirements	12V
Window	1 mm quartz
Back Focus Distance (CCD to Lens)	6.59 mm
Mounting Method	4-40 UNC
<b>Lens: Ultimate Imaging Quality under Extreme Operating Conditions</b>	
Aperture	1:1.0 in (Lowest Possible Exposure Times)
Objective Field	43 cm x 43 cm

# Appendix D - Quantum Efficiency Curves

## Kodak KAF-09000

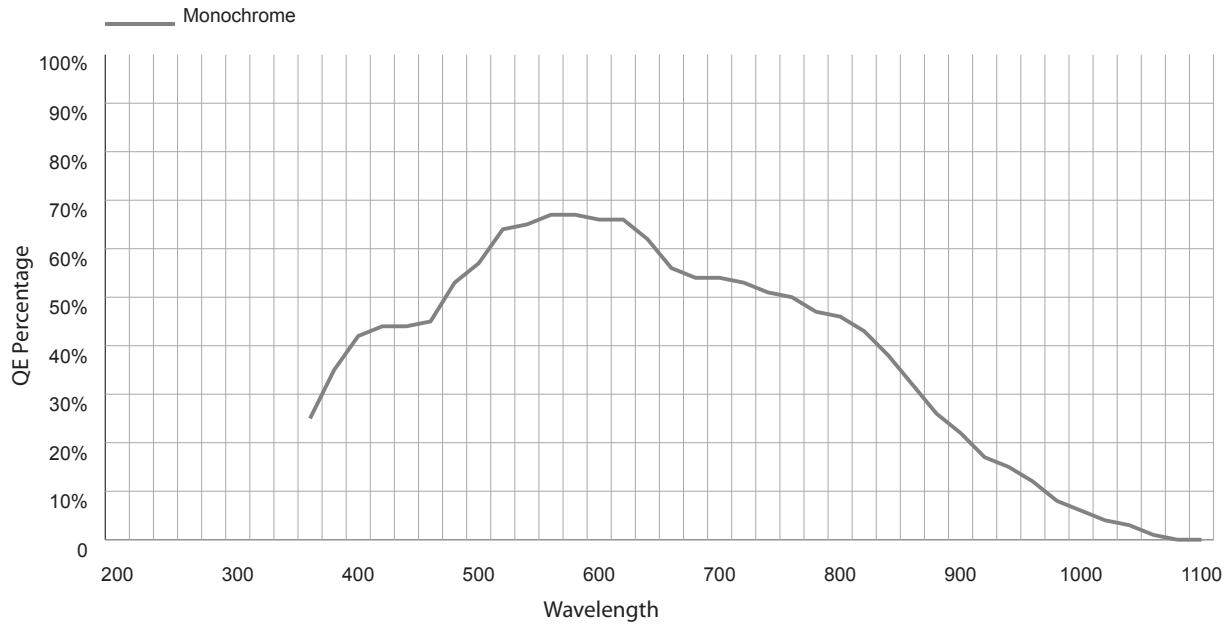

## Kodak KAF-16803

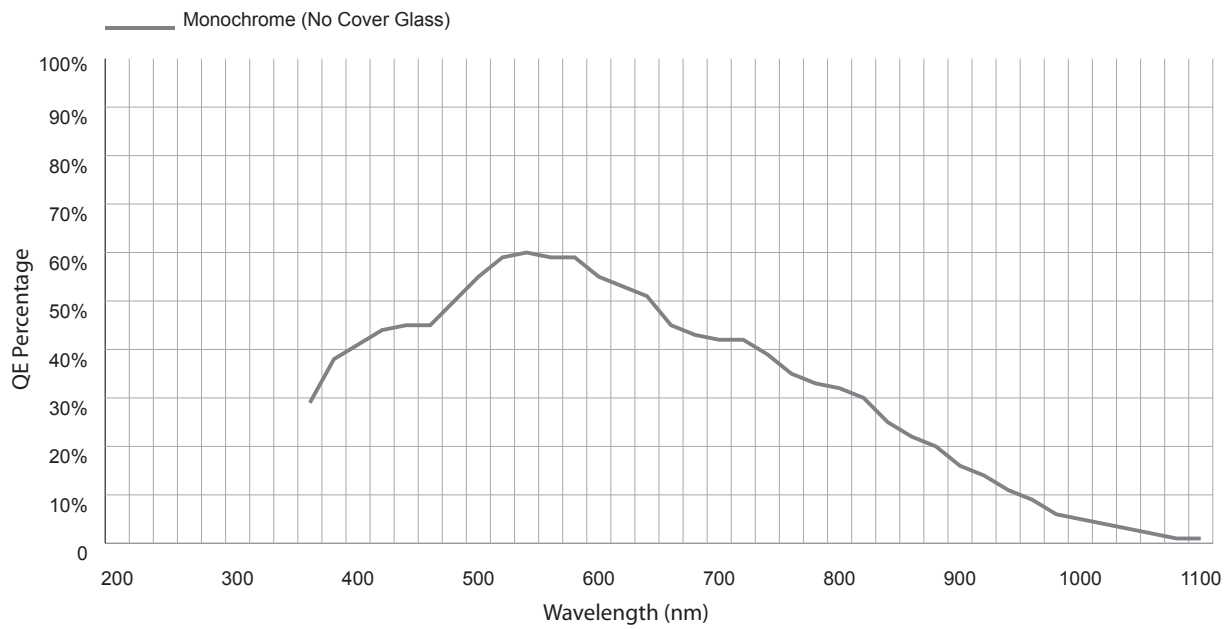

## Appendix E – Warranty for ProLine Imaging Systems

Unless otherwise noted, the standard statement of warranty described below applies to customers who purchase this System. This warranty may not apply in special circumstances in which prior arrangements have been made and separate documentation has been supplied prior to, or with, the System.

This warranty applies to ProLine Imaging Systems.

All cameras and services are FOB Lima, N.Y.. The customer is responsible for shipping and insurance to and from FLI.

The camera is warranted against defects in materials and workmanship for a period of one (1) year after delivery to the original purchaser.

The CCD array is warranted by the CCD manufacturer (Kodak) for one (1) year.

In the event of a CCD array failure or malfunction, FLI will assist in testing, replacement, shipping and required communications with the CCD manufacturer in order to facilitate a resolution of the problem.

The internal environment of the camera is warranted to remain moisture free for a period of one (1) year when used under normal conditions.

Damage arising from ESD (electrostatic discharge) events, exposure to the elements, mechanical shock, over-voltage, reverse polarity connections, or other environmental hazards is not covered under warranty.

FLI will be the sole judge of what constitutes defects vs. normal performance.

FLI application software is supplied for demonstration purposes only. The software carries no warranty of fitness for any purpose. FLI supplies the necessary information, drivers, and libraries, for users and 3rd party vendors to develop software for their specific purposes.

FLI works to maintain compatibility with many 3rd party software vendors, however FLI cannot guarantee operation with non-FLI software. FLI is not responsible for changes, upgrades, or errors in 3rd party programs.

Incidental and consequential damages resulting from the use of FLI products, malfunction or failure to perform, or lack of fitness for a particular purpose, are not the responsibility of FLI and are hereby excluded both for property damage and to the extent permitted by law, for personal injury damage.

FLI products are not authorized for use as critical components in life support or medical diagnostic applications where failure to perform could result in injury, faulty diagnosis, or other risk to patients or personnel.

FLI products are not authorized for use in robotic control systems where malfunction or failure could cause system motions hazardous to personnel.

This warranty applies to the original purchaser.

## Appendix F – Camera and Supplies Return Procedure

If you need to return a camera or supplies, please follow the instructions outlined below.

1. Obtain authorization to return the camera/products in advance by phone or email contact with FLI and:
2. If you are outside the United States, contact your Customs Authority to register the merchandise to be returned to the United States for warranty repair or refund. Use the Harmonized Code number 9801.00.1012 on your shipping documentation. The monetary value you place on the item should be stated for insurance purposes. Clearly state that the “Value is for Customs purposes ONLY.” When FLI returns the repair item to you, we will use the same monetary value.
3. For all customers, prepare a Pro Forma invoice to accompany the shipment with the following statement:
4. For Equipment not covered under warranty: “American goods returned for repair only with NO Commercial Value. Temporary return only”

OR:

- ∞ For Equipment covered under warranty: “American goods returned for Warranty Repair only with NO Commercial Value. Temporary return only”
  - ∞ C) For all customers, if you are requesting service under warranty or a return, a copy of your original receipt.
  - ∞ For you records, make a copy of these documents.
  - ∞ Prepare a large shipping label with the appropriate return address (FLI or distributor) and for shipments from outside the U.S., include the Harmonized Code number.
5. Locate the original shipping boxes in which your item(s) was packaged. These boxes are designed to protect the products.

OR:

If you do not have the original shipping boxes, obtain a rigid box that is at least 3” (7.5 cm) larger in all dimensions than the items. A smaller box will not allow appropriate cushioning. Tape the side and bottom seams to secure the box.

6. If you have the original packing materials, place the item(s) in the original plastic bag(s) and place the bagged item in the appropriate foam cutout in the proper orientation. Insert other items into their appropriate compartments.

OR:

7. If you do not have the original bag, place the item(s) in a plastic bag and seal it. Wrap the bagged item(s) with at least two layers of bubble wrap or two bubble wrap bags. Wrap other items in the same manner. Into the bottom of the box, place two inches of packing material (Styrofoam peanuts or additional bubble wrap). Place the item(s) on the bottom layer with space around each. Add additional packing material around the sides of each item(s) and on top of the item(s).
8. Write a letter that includes the following: reason the item is being returned to FLI or distributor, your complete contact information (name, phone number(s), email address, return shipping address), and if appropriate, payment method and information. 4. On top of the item(s) in the box, add the required paperwork described in step 1 and the letter described in step 4. Seal the box with packing tape. Tape the top flaps and label the box with the shipping label prepared in step 1.
9. Contact a shipper for pickup or bring it to a reliable carrier. As noted in step 1, use the appropriate value on shipping forms. FLI is not responsible for damage to any item or items when they are in the possession of a carrier.

# Return Addresses by Carrier

## **UPS and Fed Ex Returns**

---

Finger Lakes Instrumentation  
Att. Greg Terrance  
7287 West Main St.  
Lima, N.Y. 14485

## **USPS Returns**

---

Finger Lakes Instrumentation  
Att. Greg Terrance  
P.O. Box 19A  
7298 West Main St.  
Lima, N.Y. 14485